



## SK-WP Series ADDENDUM

This addendum is to the SK-WP series instruction manual.

### Installing Batteries

Battery cover lock

- To remove the battery cover, unlock the battery cover lock first.  
Be sure to lock the battery cover lock again after installing batteries.



### Automatic Power-Off Function


- To disable or enable the automatic power-off function, press the [ON/OFF] key while pressing the [RE-ZERO]/[TARE] key. The scale memorizes this setting and it will be shown by pressing both the [RE-ZERO]/[TARE] and [UNITS]/[ZERO] keys in the weighing mode.

PoFF0

Function disabled.

PoFF1

Function enabled.

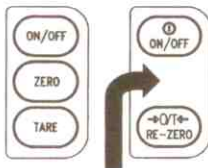

Location of the [UNITS] key for the (k)g models

### Calibration using a Calibration Weight

- ° CAL F in the instruction manual is shown as ° FULL.

- "Calibration By Gravity Compensation" is as follows.

- When CAL is displayed, press the [UNITS]/[ZERO] key.
- 97.98 The previously set value will be displayed. Press the [RE-ZERO]/[TARE] key to change the first digit after ".", ". Press [UNITS]/[ZERO] to shift the ".", and select the digit to be changed.
- 980.1 Set the new value, and press and hold the [UNITS]/[ZERO] key until End is displayed.
- After showing CAL, press [ON/OFF] to turn the power off.

### SPECIFICATIONS


Display update    Approximately 10 times per second

# AND

A&D Company, Limited

## SK-WP Scales Class III

This addendum describes about the way to enter calibration mode for the class III approved SK-WP scales.

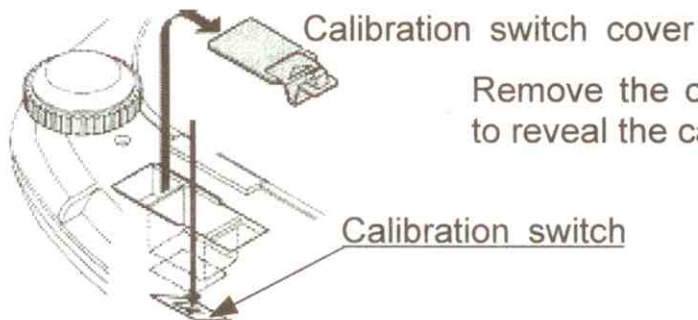

Remove the calibration switch cover to reveal the calibration switch.

## Enter Calibration Mode

Ensure that the scale is switched on and then press the calibration switch. The display will show CAL.

Note) The other procedure to calibrate the scale is the same with the standard SK-WP (not trade approved). Please refer to the attached instruction manual.